LABORATORY MINI-DRYER

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for Drying/Curing/Setting/Baking and Thermofixation Tests

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A DRYING, CURING, FIXATION APPARATUS & SETTING, BAKING AND THERMOFIXATION

Machine Designation

Having gained long term experience, RAPIDCO develop and produce the MINI-DRYER which now becomes an indispensable apparatus in laboratories to perform and assist dyeing/finishing tests and thermosoling processes.

The "RAPID" MINI-DRYER allows tests to be carried out for recipe evaluation during resin finishing, chemical tests, production for hand feeling samples, and to check degree of colour change during dyeing and finishing processes and to prepare fabric samples with the correct colour shade and structure.

The MINI-DRYER has special designs of the sample feeding device and its holding pin-frame. It is important when feeding in test samples automatically, and with a small entry to the drying chamber, that precision in temperature control should be assured, unlike conventional models which have a temperature tolerance of ± 10°C or above.

With the sample holder, the pin frame can be stretched dimensionally in all directions to meet requirements as in the pilot plant. This of course offers more advantages with knitted fabrics or other delicate samples with much elongation characteristics.

Special Features

- (1) Automatic feeding and delivery of test samples, control of dwell time also automatic.
Automatic Feeding-Time Control - Delivery (with Call signal).
- (2) Precision in Temperature Control.
 - Because of a small entry to the chamber, inside with hot-air circulation, temperature constant is maintained very stably and will not be affected when feeding in test samples.
 - Hot air applies directly to the sample through nozzles which therefore ensures precise temperature control and distribution.
Maximum working temperature 250°C (tolerance ± 1%)
- (3) Specially-designed pin frame holds all types of sample fabrics, both woven and knitted. It offers more advantages to those with much elongation characteristics and can resume the correct structure by pulling with the pin frame then results in a smooth and accurate piece.

Applications

- (1) To perform baking and thermofixation tests.
- (2) To perform and evaluate the degree of colour change and to further avoid it happens during resin finishing and/or drying processes after dyeing.
- (3) To perform dyestuff tests, auxiliary and chemical tests, and to offer similar conditions as the pilot plant to evaluate influences on fastness and product quality.

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
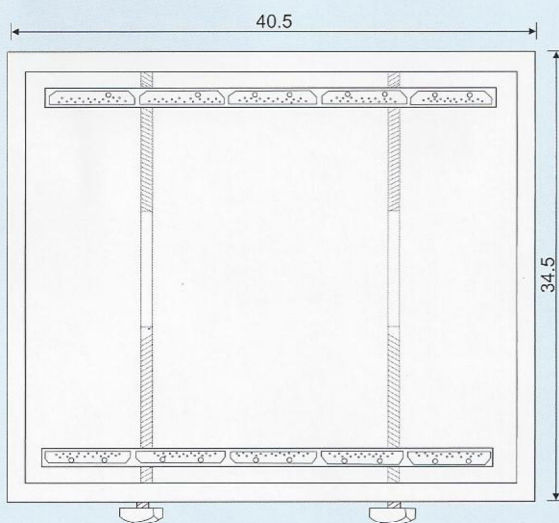
- (4) To evaluate precise resin finishing and its topping formulation.
- (5) To assist selection of chemicals and auxiliaries and to determine their concentration and quality.
- (6) To offer exactly similar conditions as for temperature, time control, and with equal elongation to prepare feeling samples to facilitate production of the pilot plant and also assist further Research & Development.
- (7) To perform water repellent, flame resistant and/or coating tests as required.
- (8) To offer a smooth and accurate test sample for higher precision obtained with computerized colour systems, to give to customers the correct fabric structure and precise and smooth colour cards.

Technical Specifications

- (1) Sample Size: 36cm \times 42cm (Maximum)
- (2) Temperature Control
 - a. Full Range Automatic Temperature Controller (Accuracy: $\pm 1\%$)
 - b. Working Temperature: 20°C to 250°C
 - c. Electric Heating: 6 kw (8 kw optional)
- (3) Dwell Time: from 10 seconds to 6 min. Preset for Automatic delivery.
- (4) Fan Motor: 3 Phase
- (5) With Damper to adjust air velocity through nozzles.
- (6) Pin Frames: Single direction latitudinal stretching 2 sides (Fig. 1)
All directions dimensional stretching to 4 sides allows 30% elongation (Fig. 2)
*Special design of pin frames available on request.
- (7) Electric connection: 380V \times 3Phase
- (8) Machine Dimensions: (76+56) \times 66 \times 80cm (L \times W \times H)

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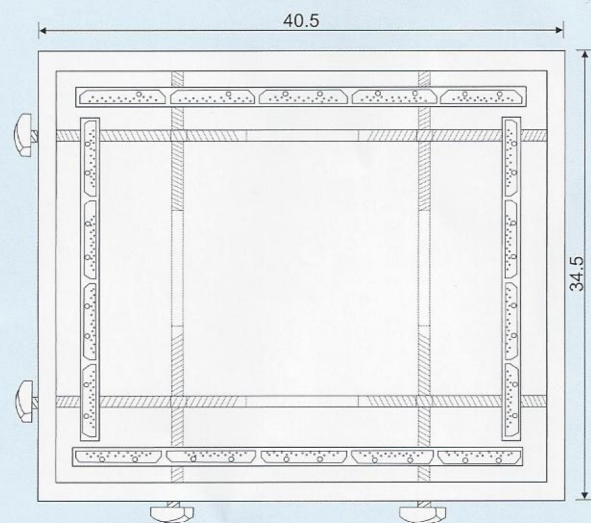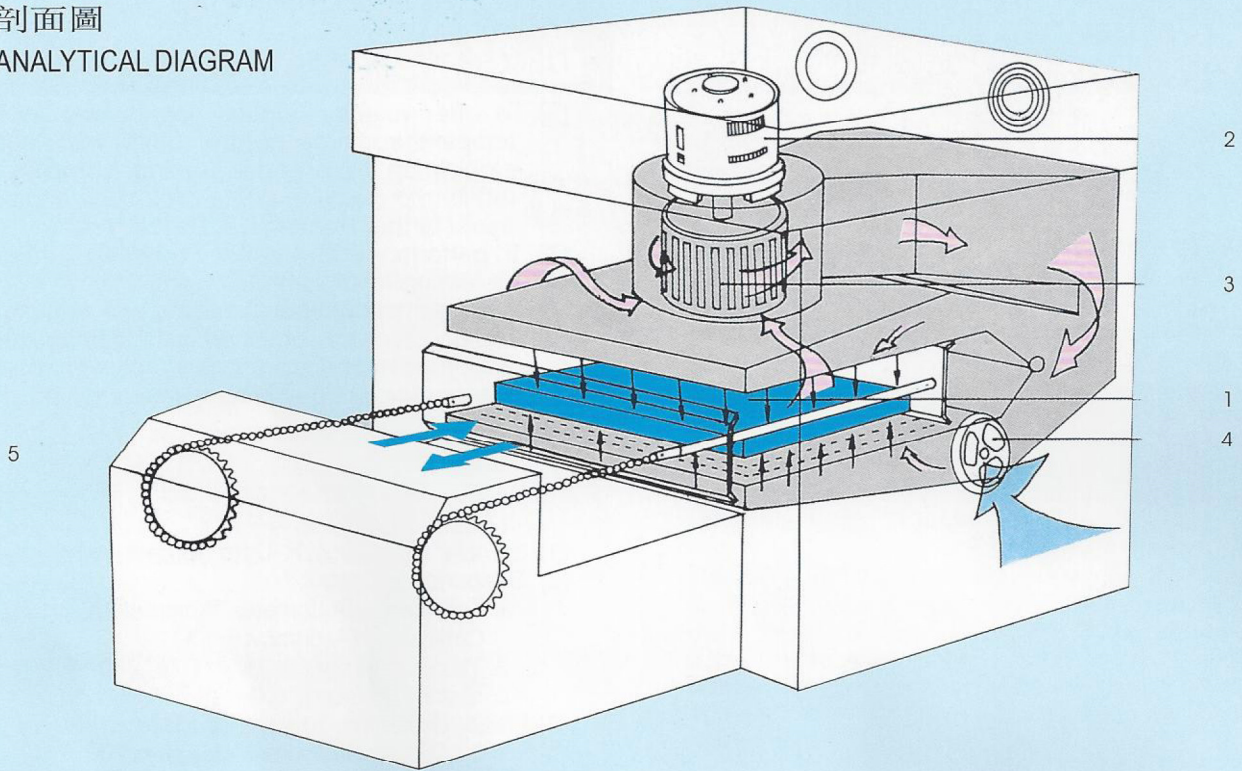
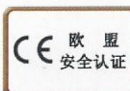
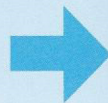
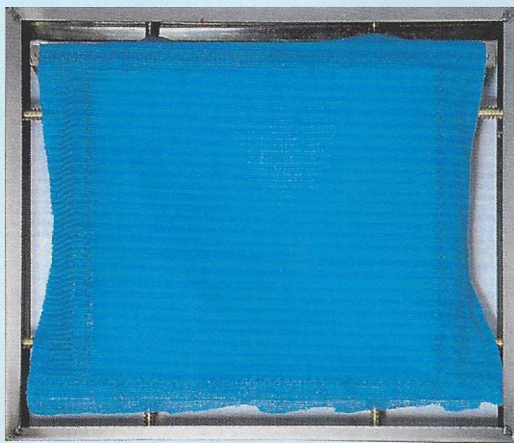
SPECIAL PIN FRAMES developed by "LABORTEX" enable precise stretching test samples to their required dimensions.

Laborteks Analitik ve Kalite Kontrol Test Cihazları San.Tic.Ltd.şti

İstoç 31.Ada No: 48/50 P.K : 34217Bağcılar / ISTANBUL

Tel : +90 212 623 00 00

Fax : +90 212 623 05 17

Gsm : +90 532 257 15 04

Web : www.trrapid.com

e-mail: apetek@laborteks.com.tr